

Frozen Cell Pellet Submission Instructions for gDNA Extraction

- Notify the WiCell Characterization lab that you will be shipping frozen cell pellets for testing (characterization@wicell.org) at least 24 hours in advance.
 - Include tracking information, if available.
 - o gDNA extraction of frozen cell pellets is available for STR and Microarray testing.
- Complete one Characterization Testing Request form for each sample and send form(s) in your shipment and/or email them to characterization@wicell.org.
- Cell pelleting protocol is available upon request.
- Frozen Cell Pellet Requirements
 - o Pellet approximately 2.0x10⁶ cells in a 1.5mL microcentrifuge tube.
 - Label sample tubes with cell line name and passage number, as you would like them to appear on the final report, and number of cells.

Frozen Cell Pellet Shipping Instructions for gDNA Extraction

- Place the 1.5mL microcentrifuge tubes within a 15mL conical tube and pack securely in a Styrofoam box.
- For international submissions, we recommend that you include a temperature monitoring system in your shipment.
- Deliveries are accepted Monday through Friday from 8:00am to 4:00pm (Please <u>DO NOT</u> ship First Overnight).
- Place Styrofoam box in a cardboard box, label, and ship OVERNIGHT or INTERNATIONAL PRIORITY to:

WiCell Characterization Lab 504 S Rosa Rd, Ste. #101 Madison, WI 53719

Shipping materials are available upon request

Please contact us at (608) 316-4145 or email characterization@wicell.org with any questions.